



# Unistat<sup>®</sup> 615w

## Ramping a 250-litre Buchi Glas Uster GLSS reactor

#### Requirement

This case study shows the remarkable power transfer capabilities of the Unistat range in using a Unistat 615w to heat and cool a 250-litre Buchi Glas Uster GLSS reactor.

#### Method

The Unistat was connected to the reactor using two 2-metre insulated metal hoses. The reactor was filled with 200 litre of Ethanol.

#### Results

The closeness of control can be seen as the Unistat ramps the process temperature between 20 °C to 0 °C and back at differing preprogrammed ramp rates.

### Setup details

Temperature range: Cooling power: Heating power: Hoses: HTF: Reactor:

-60...200 °C 9.5 kW @ 200...0 °C 8.0 kW @ -20 °C 4.8 kW @ -40 °C 1.2 kW @ -60 °C 12 kW M38x1,5; 2x2 m DW-Therm Buchi Glas Uster CR252 250-litre glass-lined (enameled) steel reactor 200 litre Ethanol 1: 90 rpm process





